




LyfoCults[®] Plus

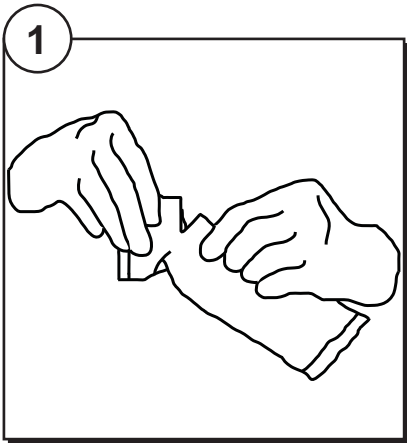


1
Tear open foil pouch at notch and remove LyfoCults Plus device

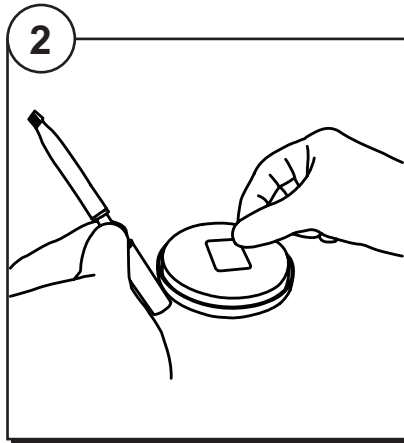


2
Remove label and place on culture media.

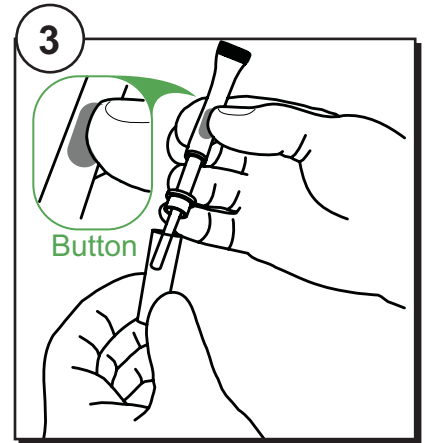


3
Unscrew applicator from vial. Holding applicator at a slight angle to vial, activate ampoule by squeezing button firmly.

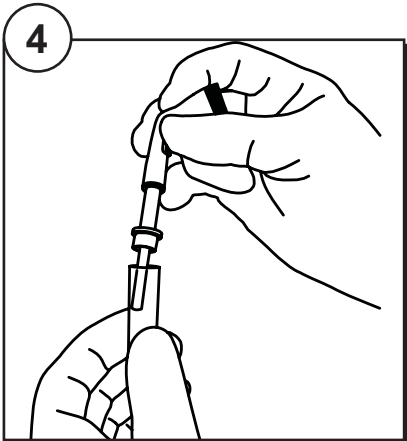


4
With the applicator at a slight angle to the vial, squeeze the ampoule dispensing the hydration fluid into vial.

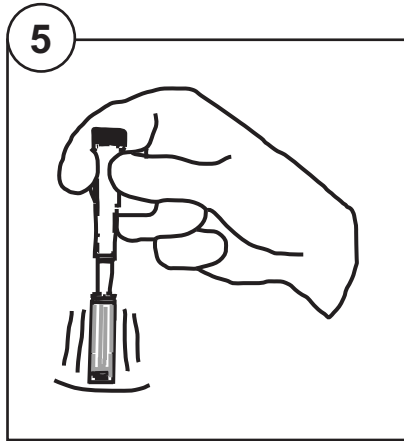


5
Re-insert applicator to vial and gently mix from side-to-side

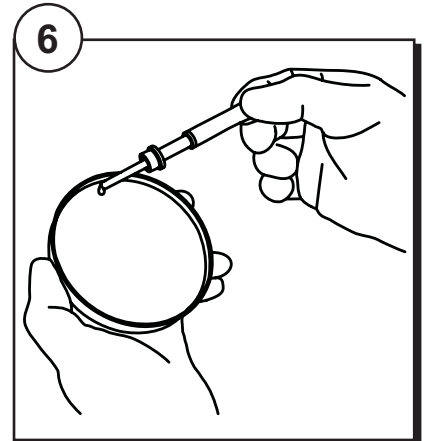

6
Using the ampoule as a transfer pipette, place a drop of resulting suspension on culture media.

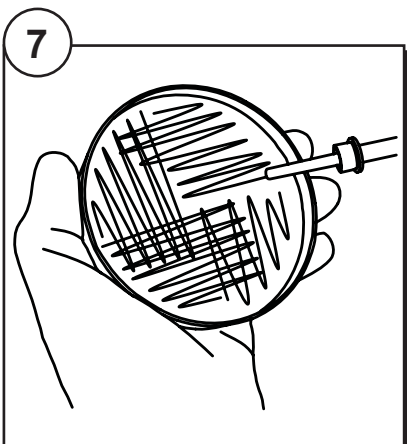

7
Streak the media for isolation and incubate in appropriate atmosphere

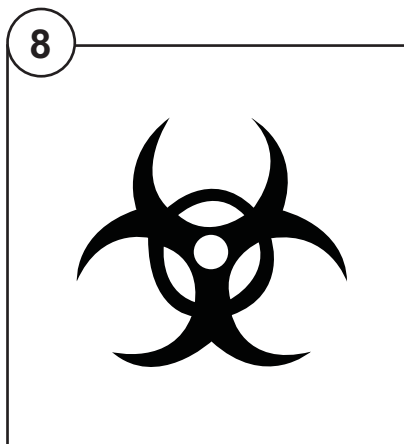

8
Discard used LyfoCults Plus appropriate biohazard receptacle

To purchase LyfoCults[®] Plus
or for further information,
please contact us

Quality Technologies, LLC

Tel: 805-498-4870

Toll Free: 800-222-0728

Fax: 805-375-7314

Web: www.quality-technologies.net